Roxy M.™

Repeating Pipettor

Catalog No. F37310-0000

Instructions for use

1. Pipettor overview

1 Dispensing

By fully pressing the dispensing button down until it stops, the adjusted volume is dispensed.

2 Adjusting wheel

By moving the adjusting wheel into the positions 1 to 5, the desired dispensing volume is set in compliance with Table 1.


3 Filling slide

The inserted tip is filled by pulling the slide upward.

4 Stopping lever

The stopping lever is used to secure the inserted tip.

5 Replaceable dispenser tip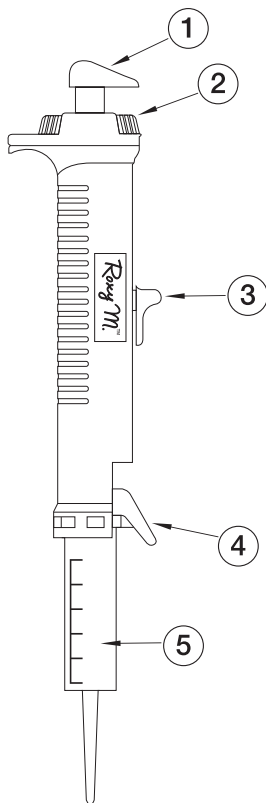

2. Selecting the volume

- Use Table 1 to determine setting for the desired dispensing volume.
- The position, to which the adjusting wheel (2) has to be set, is given in the topmost line.
- In the left-hand column, identify the tip size that will be inserted.
- Read the dispensing volume from the chart.
- For the number of dispensing doses, please refer to the bottom most line in Table 1.

Table 1


Tip Size	Adjusting Wheel Setting				
	1	2	3	4	5
0.05 ml	1 µl	2 µl	3 µl	4 µl	5 µl
0.5 ml	10 µl	20 µl	30 µl	40 µl	50 µl
1.25 ml	25 µl	50 µl	75 µl	100 µl	125 µl
2.5 ml	50 µl	100 µl	150 µl	200 µl	250 µl
5.0 ml	100 µl	200 µl	300 µl	400 µl	500 µl
12.5 ml	250 µl	500 µl	750 µl	1000 µl	1250 µl
25.0 ml	500 µl	1000 µl	1500 µl	2000 µl	2500 µl
50.0 ml	1000 µl	2000 µl	3000 µl	4000 µl	5000 µl
Steps	48	23	15	11	8

3. Inserting dispenser tip (see notes to right)

- Push the filling slide (3) down to the stop.
- Lift the stopping lever (4) upward.
- Align the neck collar of the tip with the groove in the pipettor. Exert a slight pressure to insert tip (5), which has to lock into place.
- Push down on the stopping lever (4), to lock the tip in place.
- Use the adjusting wheel (2) to set the desired dispensing volume.

Notes for step 3:

- Be sure to select the volume setting before inserting a tip as adjustment wheel setting #5 may not be accessible otherwise.
- The 50 ml tip can be inserted/locked into place in the dispenser by means of the enclosed adapter; the 25 ml tip adapter is sold separately. For this purpose, place the adapter on the tip, and fasten by turning (bayonet socket). After use of the tip, the adapter can be removed again and can be reused as often as required.
- If desired, the 0.5 ml, 1.25 ml, and 2.5 ml tips can be used with yellow 100 µl pipetting tips affixed to the ends. This will help reduce the force of dispensing to avoid backsplashes that can occur when pipetting small volumes. The 0.05 ml tip can be modified in the same way, but only by affixing a 10 µl pipetting tip.

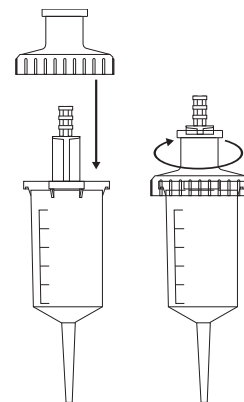4. Filling the dispenser tip (see notes to right)

- Immerse the tip into the liquid.
- Fill the tip by slowly raising the filling slide (3) upwards.
- Remove the tip from the liquid. Proceed to the dispensing steps.

Notes for step 4:

- Raise the filling slide lever slowly to the upper stop to avoid formation of air bubbles.
- An air bubble which might be in the upper part of the cylinder will have no influence on the accuracy as the last dispensing step will not be carried out due to the residual stroke blockage.
- The tips need not be filled completely.

5. Dispensing

(see notes to right)

- Wipe the tip with a lint-free tissue.
- Place the opening of the tip against the wall of the vessel.
- Dispense the liquid by fully pressing down the dispensing button until it stops, then allow it to return to its original position.
- Repeat as desired.

Notes for step 5:

- To ensure accurate dispensing, the first dispensing step must be discarded.
- The tip must be changed when the reagent is changed.
- Discard the residual liquid by slowly pushing the filling slide (3) down until it stops.

6. Removing the dispenser tip (see notes to right)

- If tip is not already empty, empty it by sliding the filling slide (3) down until it stops.
- Lift the stopping lever (4).
- Remove the tip from the dispenser and discard in accordance with local state and federal regulations.

Notes for step 6:

- When the 25 ml and 50 ml tips are used, the reusable adapters can be removed from the tip, by turning counter-clockwise, after removal from the dispenser.
- Avoid touching the outlet opening when removing the tip (beware of potentially aggressive solutions).
- When using viscous solutions, the tightness of the tip is not guaranteed after several pipettings.

7. Maintenance (see notes to right)

- The dispenser requires no maintenance or calibration. All internal settings for accurate and precise dispensing have been made in the factory. If you encounter problems, please see section 8 on troubleshooting.

Notes for step 7:

- The dispenser must not be disassembled. Such action will void the warranty.

8. Troubleshooting

Problem	Reason	Remedy
Tip cannot be inserted	Filling slide has not been lowered to the stop	Lower filling slide until it stops
Stopping lever cannot be pushed down	Tip has not been placed properly	Remove tip, align tip with groove and push tip into place
Filling slide cannot be moved upwards	Stopping lever has not been pushed down completely	Push down on the stopping lever until you hear it click into place
Dosing volume is uncertain	Adjusting wheel has not locked into place	Lock adjusting wheel into place

9. Technical Data

Weight: approx. 100 g

Accuracy and precision: refer to Table 2

Liquid: double distilled water

Reference temperature: 20° to 25 °C,
± 0.5° constant

Measurement: in accordance with German standard EN ISO 8655

Subject to technical modifications

Made in Germany

Table 2

Adjustment on dispenser	Dispensing Volumes					Systematic Error	Random Error
	1	2	3	4	5		
Number of steps	48	23	15	11	8		
0.05 ml	1 µl	2 µl	3 µl	4 µl	5 µl	± 6	≤ 9
0.5 ml	10 µl	20 µl	30 µl	40 µl	50 µl	± 1.0	≤ 1.5
1.25 ml	25 µl	50 µl	75 µl	100 µl	125 µl	± 0.9	≤ 0.9
2.5 ml	50 µl	100 µl	150 µl	200 µl	250 µl	± 0.7	≤ 0.9
5.0 ml	100 µl	200 µl	300 µl	400 µl	500 µl	± 0.3	≤ 0.8
12.5 ml	250 µl	500 µl	750 µl	1000 µl	1250 µl	± 0.3	≤ 0.7
25.0 ml	500 µl	1000 µl	1500 µl	2000 µl	2500 µl	± 0.5	≤ 0.6
50.0 ml	1000 µl	2000 µl	3000 µl	4000 µl	5000 µl	± 0.5	≤ 0.6

The data for tips 0.5, 1.25 and 2.5 ml are based on a measurement carried out with a 100 µl pipetting tip. The measurement for tip 0.05 ml was performed with a 10 µl pipetting tip. Tips and dispenser have received the certificate of conformity with the symbol below in accordance with the Weights and Measures Regulations as amended on 12th August 88.

10. Limited Warranty

Bel-Art Products Roxy M[®] manual repeating pipettors are warranted against manufacturing defects in materials and workmanship for a period of one year from the date of purchase by the original consumer or other original end user. The exclusive remedy for any and all warranties and claims related to any product of Bel-Art Products is the return of the product for replacement or repair, at Bel-Art Products' discretion except that such remedy may be further limited as outlined on www.belart.com/warranty. To activate your warranty, please complete the warranty registration within 10 days of receipt of product at www.belart.com/warranty.

Ordering Information

Description	Catalog No.		Unit of Measure	
Roxy M [®] Repeating Pipettor	F37310-0000		Each	
Roxy M [®] Tips	Standard	Sterile	Standard	Sterile
Roxy M [®] Tips, 0.05ml	F37300-0000	F37300-0017	100/Pack	100/Pack
Roxy M [®] Tips, 0.5ml	F37300-0001	F37300-0010	100/Pack	100/Pack
Roxy M [®] Tips, 1.25ml	F37300-0002	F37300-0011	100/Pack	100/Pack
Roxy M [®] Tips, 2.5ml	F37300-0003	F37300-0012	100/Pack	100/Pack
Roxy M [®] Tips, 5.0ml	F37300-0004	F37300-0013	100/Pack	100/Pack
Roxy M [®] Tips, 12.5ml	F37300-0005	F37300-0014	100/Pack	100/Pack
Roxy M [®] Tips, 25ml	F37300-0008	F37300-0015	25/Pack	25/Pack
Roxy M [®] Tips, 50ml	F37300-0009	F37300-0016	25/Pack	25/Pack
Roxy M [®] Tips Assortment	F37300-0100	-	45/Pack	-
Roxy M [®] 25ml Tip Adapter - Red	F37300-1025	-	3/Pack	-
Roxy M [®] 50ml Tip Adapter - Gray	F37300-1050	-	3/Pack	-

661 Route 23 South, Wayne, NJ 07470

TEL: 973-694-0500 • FAX: 973-694-7199 • www.belart.com

Roxy M[®] is a Trademark of Bel-Art Products.

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

© 2020. All Rights Reserved. Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.

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© 2020. All Rights Reserved. Bel-Art® - SP Scienceware® is a division of SP Industries, Inc.
© 2020. All Rights Reserved. Bel-Art[®] - SP Scienceware[®] is a division of SP Industries, Inc.