Instrucciones

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BATTERY INSTALLATION

1. Use a suitable size coin to open the battery cover.
2. Install the battery with correct polarity positioning.
3. Close the battery cover.

Place the probe in the bottle, place bottle with probe in test condition, record your reading when the reading becomes stable.

POWER ON/OFF

Press [ON/OFF] button to switch on or off the unit.

OPERATION

1. °C / °F Exchange
Press [°C/°F] button to change temperature unit.
2. Maximum/minimum reading memory
a. Press [Max/Min] button to display maximum measured value.
b. Press [Max/Min] button once more to display minimum measured value.
c. Press [Max/Min] button again to return to normal display.
d. Press and hold [Max/Min] button for about 3 seconds then release to reset memory reading to current value.

If the display becomes dim then change the battery with a same type new battery.

Battery Handling/Packaging

- Remove batteries from equipment immediately after they fail to operate the equipment.
- If equipment uses two or more batteries, always replace batteries in complete sets.
- Do not attempt to recharge alkaline batteries, they are not rechargeable.
- Do not handle hot or warm batteries.
- Do not heat, incinerate, crush, puncture or mutilate batteries.
- Wear personal protective equipment if batteries show signs of leakage, bulging, swelling or deformation.

Warning: If potassium hydroxide electrolyte from an alkaline battery comes in contact with the skin, do not try to neutralize the electrolyte with vinegar or any other acidic solutions. Neutralization may trap electrolyte on the skin. Flush the affected skin area with copious amounts of water. If the battery electrolyte gets into your eyes, it can cause severe damage and/or blindness.

To register this product for a one-year limited warranty, visit www.belart.com/warranty

H-B Instrument – SP Scienceware
Triple Accredited/Registered
ISO 9001:2008 Registered Manufacturer
ISO 14001:2004 Registered Manufacturer
ISO/IEC 17025:2005 Accredited Laboratory

For more information
973-694-0500 (p) • 973-694-7199 (f)
info@belart.com • www.belart.com/hb

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