Instrucciones

INSTALACIÓN

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Instructions

DURAC® Single Channel Electronic Timer

INSTALLATION

Open the battery compartment by sliding the battery compartment cover

Install 1 AAA batteries following the correct polarity

Close the battery compartment by sliding the battery cover on

The display will show "00:00"

OPERATION

Count Down Function

Press the [MIN] button to set the minutes to count down

Press the [SEC] button to set seconds to count down

Press the [START/STOP] to start the timer counting down

The alarm will go off when the display counts down to 00:00

Press the [START/STOP] to stop counting - the remaining time will be displayed

Press the [START] to continuing the counting

Press the [MIN] and [SEC] buttons simultaneously to reset the display value to "00:00"

Count Up Function

Press the [START/STOP] to start the timer counting up

Press the [START/STOP] to stop counting - the elapsed time will be displayed

Press the [START/STOP] to continuing the counting

When display counts up to maximum range the timer will stop

Battery Handling/Packaging

- Remove batteries from equipment immediately after they fail to operate the equipment.
- If equipment uses two or more batteries, always replace batteries in complete sets.
- Do not attempt to recharge alkaline batteries, they are not rechargeable.
- Do not handle hot or warm batteries.
- Do not heat, incinerate, crush, puncture or mutilate batteries.
- Wear personal protective equipment if batteries show signs of leakage, bulging, swelling or deformity.

Warning: If potassium hydroxide electrolyte from an alkaline battery comes in contact with the skin, do not try to neutralize the electrolyte with vinegar or any other acidic solutions. Neutralization may trap electrolyte on the skin. Flush the affected skin area with copious amounts of water. If the battery electrolyte gets into your eyes, it can cause severe damage and/or blindness.

To register this product for a one year limited warranty, visit www.belart.com/warranty

H-B Instrument – SP Scienceware
Triple Accredited/Registered
ISO 9001:2008 Registered Manufacturer
ISO 14001:2004 Registered Manufacturer
ISO/IEC 17025:2005 Accredited Laboratory

For more information
973-694-0500 (p) • 973-694-7199 (f)
info@belart.com • www.belart.com/hb

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