


# INSTRUCTIONS



## Vacuum-Pressure Indicator

Catalog No. H42042-0000

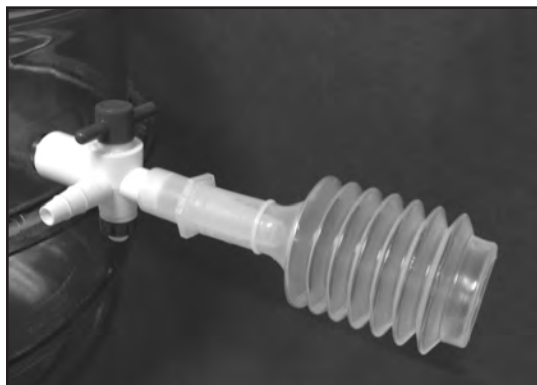

Position to pull vacuum



Position when closed

The Vacuum-Pressure Indicator, when connected to a desiccator, clearly shows when the desiccator is under vacuum.

1. Remove red plug from end of stopcock.
2. Place the Vacuum-Pressure indicator onto the tapered smooth outlet where plug was removed from.
3. When vacuum is applied to the stopcock the bellows will collapse. When the stopcock is closed, and the vacuum source removed, the bellows remains collapsed as long as there is a vacuum in the desiccator.
4. If the desiccator loses vacuum, the bellows will expand to alert the operator.

## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

942042001

05/06


Scienceware®, is a reg. TM of Bel-Art Products  
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

## OTHER FINE SCIENCEWARE® PRODUCTS FOR YOUR LABORATORY

### *Demand More From Your Desiccator!*



Secador™ 1.0  
Volume .75 cubic feet

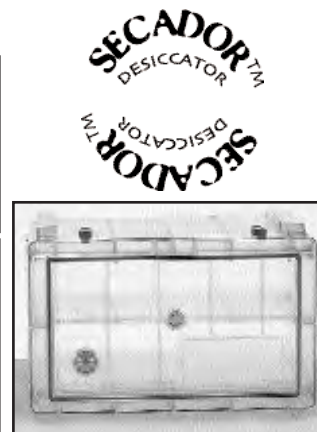

Secador™ 4.0 Horizontal & Vertical  
Volume 2.0 cubic feet

### Select the Secador™ Desiccator Cabinet that suits your lab!

- Patented sealed construction – Airtight, dust and moisture free storage
- Heavy duty copolyester composition builds rugged cabinet with long service life
- Imbedded UV protection blocks 99% of light from cabinet and contents
- Superior Chemical Resistance - Prevent staining and crazing
- Built in Hygrometer for easy RH monitoring
- Stackable – Saves valuable bench space
- Use with any Scienceware® desiccant

Description	Catalog Number
Secador™ 1.0 Desiccator Cabinet	42071-0000
Secador™ 4.0 Desiccator Horizontal Cabinet	42074-0000
Secador™ 4.0 Desiccator Vertical Cabinet	42074-1000



## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com