



Dry Keeper™ Plus Desiccator Cabinet

Catalog Number H42061-0000

To ensure that your items are being placed into a dry environment, we suggest that your desiccator be allowed to operate while empty for 24 hours.

Warning: *Do not use this cabinet to store volatile or flammable materials*



Operation

1. Place the cabinet on a level surface and plug into a 115V outlet to start the continuous automatic drying cycles. During operation, there will be normal sounds from the timer and shutters inside the drying module.
2. Adjust shelf height to accommodate items and place items on shelves
3. For maximum long term drying performance, minimize the amount of time that the door remains open and do not disconnect the cabinet from the electrical source.
4. To clean the cabinet, use a soft cloth. Do not use abrasive cleansers which will leave scratches. We recommend the use of Aquet™ Detergent, Catalog No. F17094-0030 or Cleanware™ Plastic Laboratory Cleaner Catalog No. F17084-0000.

BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

N O T I C E !

Desiccator Cabinets

H42056-0001, H42056-1003, H42058-0001, H42058-0003, H42061-0000

The hygrometer supplied with this desiccator cabinet is not a precision instrument. It registers high or low humidity, but it is not an accurate indicator of the percent of relative humidity shown on the dial.

If the hygrometer reading becomes lower after the unit has been plugged in for one complete cycle (5 hours or more), the drying mechanism is working. If a more precise measure of the actual relative humidity in the cabinet is required, a portable hygrometer may be placed in the cabinet.

661 Route 23 South, Wayne, NJ 07470

TEL: 1-800-423-5278 • FAX: 973-694-7199 • www.belart.com



942056001 – 05/15

N O T I C E !

Desiccator Cabinets

H42056-0001, H42056-1003, H42058-0001, H42058-0003, H42061-0000

The hygrometer supplied with this desiccator cabinet is not a precision instrument. It registers high or low humidity, but it is not an accurate indicator of the percent of relative humidity shown on the dial.

If the hygrometer reading becomes lower after the unit has been plugged in for one complete cycle (5 hours or more), the drying mechanism is working. If a more precise measure of the actual relative humidity in the cabinet is required, a portable hygrometer may be placed in the cabinet.

661 Route 23 South, Wayne, NJ 07470

TEL: 1-800-423-5278 • FAX: 973-694-7199 • www.belart.com



942056001 – 05/15

N O T I C E !

Desiccator Cabinets

H42056-0001, H42056-1003, H42058-0001, H42058-0003, H42061-0000

The hygrometer supplied with this desiccator cabinet is not a precision instrument. It registers high or low humidity, but it is not an accurate indicator of the percent of relative humidity shown on the dial.

If the hygrometer reading becomes lower after the unit has been plugged in for one complete cycle (5 hours or more), the drying mechanism is working. If a more precise measure of the actual relative humidity in the cabinet is required, a portable hygrometer may be placed in the cabinet.

661 Route 23 South, Wayne, NJ 07470

TEL: 1-800-423-5278 • FAX: 973-694-7199 • www.belart.com



942056001 – 05/15